

Ringbio


How to use the incubator

① ----- ② ----- ③ ----- ④ ----- ⑤

① turn on your incubator, your screen will like this.



② set the temperature, for example, 40°C. Short press the “Prog”, then move to the temp position by pressing “left” or “right”, then set wanted temp by pressing “up” or “down”.

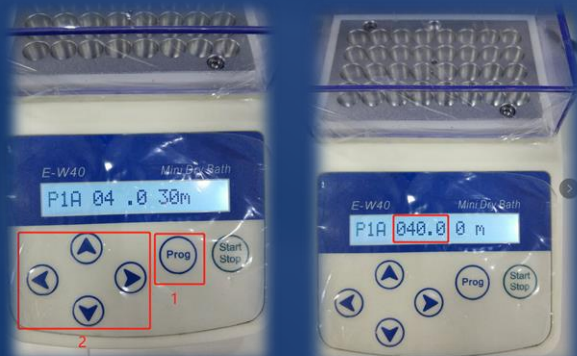

③ set the time, for example, 3min. Move to the time position by clicking “left” or “right”, then set wanted time by pressing “up” or “down”.



④ **long press** the “Prog” to finish the setting, then wait the incubator reach up to your setting temperature, finally, the screen will like this.



⑤ put your well(reagents) or test strip on the incubator, short press “Start” to run the program.

Ring Biotechnology Co., Ltd

E-mail: export@ringbio.com diego@nbgen.com Web: www.ringbio.com

Add: Building 3, Zhongtongtai TechnoPark, No. 11, Kechuang 14th St, Beijing 100176, CHINA


Tel: +86-10-56267496 Technical Support & Service: +86-18600362934

Ringbio


How to conduct a standard test




1 Bring all test samples to room temperature; number them to keep record.

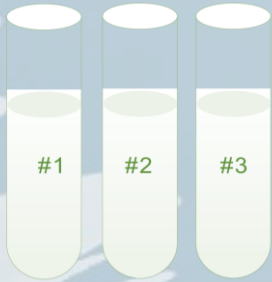


2 Take test kit according to your sample number and also number the kit wells to keep record and consistency.

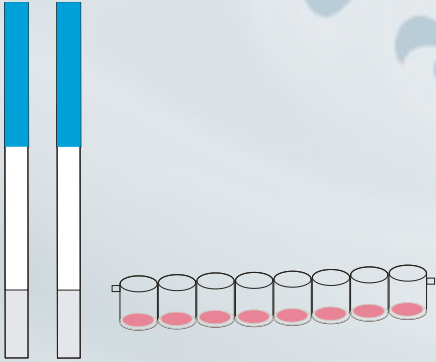


3 Take 200ul sample into the wells using pipet. You can also then put the well into the well holder to avoid sample spill.

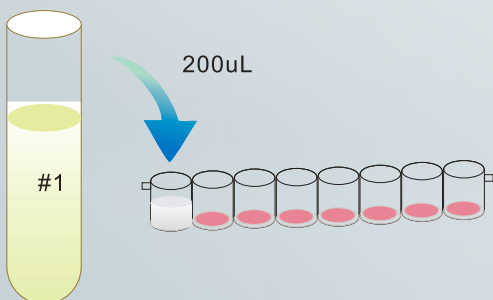

4 Absorb up and down for 5 times to mix sample with reagent completely. Start the timer when the mixture is pink. Incubate at 40 °C for 3 min.

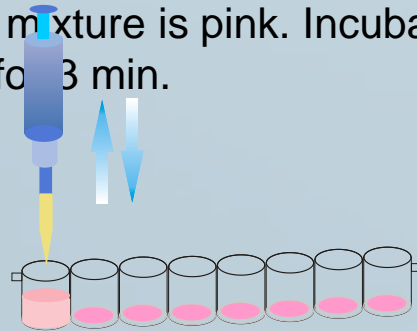

5 Insert the "Immersed" end of the strip into the mixture; Incubate for 3 min at 40 °C again.

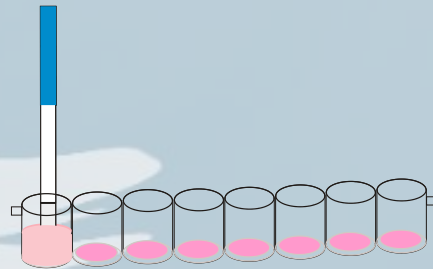

6 Take out the strip; judge the result according to kit instruction.

